



Food Packaging Residual Solvents on MEGA-VOC 1 & MEGA-VOC 2 dual system

MEGA-VOC 1 & 2 kit available for single injector port,
 dual detector GC system.

Contact us for more information.

Conditions:

Injection: Split 250°C, 0.2µL with micro-syringe, 150mL/min Split Flow.

Detector: FID 250°C.

Oven Temp.: 35°C (0.4 min), 35°C/min, 200°C.

Carrier Gas: Helium, 60 kPa.

Peaks:

1. Ethanol
2. Acetone
3. 2-propanol
4. 1-propanol
5. Methyl Ethyl Ketone
6. Ethyl acetate
7. Cyclohexane
8. Isobutanol
9. Isopropyl acetate
10. 1-butanol
11. Methoxy propanol
12. Propyl acetate
13. Isobutyl methyl ketone
14. Ethoxy propanol
15. Toluene
16. Butyl acetate

